BRESSER[®]




USB Hand Microscope LCD

Art. No. 88-54200

- Ⓛ DE Bedienungsanleitung
- Ⓛ GB Operating Instructions
- Ⓛ FR Mode d'emploi
- Ⓛ NL Handleiding
- Ⓛ IT Istruzioni per l'uso
- Ⓛ ES Instrucciones de uso
- Ⓛ PT Manual de utilização

DE	Bedienungsanleitung	4
GB	Operating Instructions	18
FR	Mode d'emploi	30
NL	Handleiding	44
IT	Istruzioni per l'uso	58
ES	Instrucciones de uso	72
PT	Manual de utilização	86

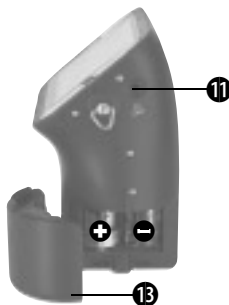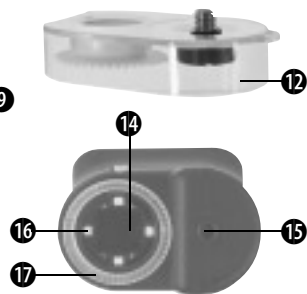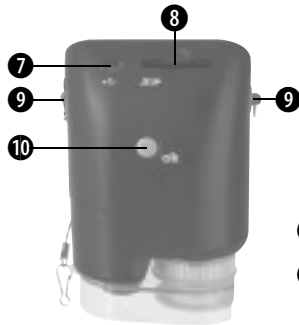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RISK to your child.

 This device contains electronic components that are powered by either a mains connection or batteries. Never leave a child unsupervised with this device. The device should only be used as per these instructions otherwise there is a serious RISK of ELECTRICAL SHOCK.

Batteries should be kept out of children's reach. When inserting batteries please ensure the polarity is correct. Leaking or damaged batteries can cause injury if they come into contact with the skin. If you need to handle such batteries please wear suitable safety gloves.

Children should only use this device under supervision. Keep packaging materials (plastic bags, rubber bands, etc.) away from children. There is a risk of SUFFOCATION.

FIRE-/ DANGER OF EXPLOSION!

 Do not expose the device to high temperatures. Use only the mains adapter supplied or those battery types recommended. Never short circuit the device or batteries or throw into a fire. Exposure to high temperatures or misuse of the device can lead to short circuits, fire or even explosion!

RISK of material damage.

 Never take the device apart. Please consult your dealer if there are any defects. The dealer will contact our service centre and send the device in for repair if needed.

Do not subject the device to temperatures exceeding 60 C.

TIPS on cleaning

Remove the device from its energy source before cleaning (remove plug from socket / remove batteries)

Clean the exterior of device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components.

Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging. Batteries should be removed from the device if it is not going to be used for a long period of time.

DISPOSAL

Dispose of the packaging material/s as legally required. Consult the local authority on the matter if necessary.

Do not dispose of electrical equipment in your ordinary refuse.

The European guideline 2002/96/EU on Electronic and Electrical Equipment Waste and relevant laws applying to it require such used equipment to be separately collected and recycled in an environment-friendly manner.

Empty batteries and accumulators must be disposed of separately. Information on disposing of all such equipment made after 01 June 2006 can be obtained from your local authority.

EEC conformity explanation

Meade Instruments GmbH & Co. KG has formed a “Conformity Declaration” in line with the applicable guidelines. This declaration is available for review.

These are the parts of your microscope

1. Optical Zooming switch
2. Power LED
3. POWER & ZOOM IN & MENU button
4. LCD monitor
5. FUNCTION (Right) button
6. PLAYBACK (left) button
7. USB Port
8. SD Card Socket
9. Strap Eyelet
10. SHUTTER/OK button
11. TV OUT (optional) port
12. View case (optional)
13. Battery Compartment Cover
14. Lens
15. Tripod /Screw Socket
16. LED lights
17. Magnification Revolving Switch
18. Neck strap
19. Wrist strap
20. Video Cable (optional)
21. Software CD and USB Cable

1. Install Batteries

1. Rotate the screw. Take off the View case.
2. Slide and open the battery compartment cover.
3. Put in the batteries in the correct direction. + and - terminals has shown..
4. Slide and close the battery compartment cover.

2. Driver and Software Installation

2.1 Installing the Drivers

Attention!

Do not connect the „USB Hand Microscope LCD“ to your PC (computer) yet. In order to work with the „USB Hand Microscope LCD“ for the long term, we recommend that you always connect it to the same USB port. In some cases, your computer may not detect the device. If this occurs, you can typically fix the issue by re-installing the drivers

(from the CD). If this does not work, please see the chapter on troubleshooting.

Insert the supplied CD-ROM into your CD/DVD drive. The CD menu opens automatically. If the menu does not open, select the CD/DVD drive in Windows Explorer. There, start the file „Menu.exe“ with a double click.

Please make sure that the „USB Hand Microscope LCD“ is not connected to your PC.

Take the following steps:

1. Save the current data and close any open programs.
2. Click on „Treiber Installieren“.
3. The „Setup Status“ window now appears. The program files are copied to your PC. This process can take a few minutes.
4. After the window disappears, it can take a few moments before the „InstallShield Wizard Complete“

window appears. Click on „Finish“.

5. The driver is installed now.

2.2 Installing the Image Editing Software

Note: If you have already installed an image editing program, you do not need to go through with this Point. In this way, only the driver for the „USB Hand Microscope LCD“ is installed. If you are not sure, please continue with the following steps.

1. Insert the supplied CD-ROM into your CD/DVD drive. The CD menu opens automatically. If the menu does not open, select the CD/DVD drive in Windows Explorer. There, start the file „Menu.exe“ with a double click.

2. Click on "Software Installieren"

3. A selection menu appears. Here, you can select the desired language. Confirm with „OK“.

4. Click „Continue“ in the „Welcome“ screen.

5. The License Agreement appears. Confirm with „Yes“.

6. The next two windows will ask you for the „Destination Directory“ and the „Program Folder.“ Confirm this with „Continue“.

6. The „Setup Status“ window appears. The program files are copied to your PC. This process can take a few minutes.

8. The „InstallShield Wizard Complete“ window appears. Click on „Finish“. The image editing software is now installed.

2.3 Troubleshooting: the device isn't recognised by the computer (USB)

1. This occurs in rare cases with some OEM Windows Vista or XP versions. The device is not defective. Updating the notebook or PC USB driver usually suffices.

2. If not, proceed as below. (The maker cannot accept any liability whatsoever in these cases¹).

Delete ghost devices. These are devices not actually currently connected to your computer. The reason for this is that Windows generates a system settings entry for each USB device (e.g. USB Stick) on each port. It then seeks that/those device/s each time it's started. This slows down USB device recognition and Windows booting and causes non-recognition of USB ports.

Cleaning up the system environment and deleting devices that are no longer in use helps here. To do so right click WORKPLACE (COMPUTER in Vista) and open the PROPERTIES menu. Then click the ADVANCED (ADVANCED SYSTEM SETTINGS in Vista) and then ENVIRONMENT VARIABLES. Choose the NEW option in ENVIRONMENT VARIABLES. Enter the following under VARIABLE NAME.

devmgr_show_nonpresent_devices
Under VARIABLE VALUE enter „1“. Confirm your input with OK and then

reboot the computer.

Change to the device manager. In the VIEW menu activate the SHOW DEVICES OPTION. The ghost devices previously not shown will now be displayed in light grey. Check the various categories including USB, memory size, etc. Only delete entries for devices you no longer use.¹

¹Source: Microsoft Knowledge Base <http://support.microsoft.com/kb/315539/de>

3. Using the Microscope

3.1 Insert an SD card:

1. Turn off the microscope.
2. Insert the card in the correct direction as shown. Push the card into the slot and be locked.
3. To remove the card, push it in, then release it. When the card is partially ejected, pull it out.

Note: If the “MEMORY FULL” warning message appears on the

LCD monitor, it means the SD card is full. Please transfer all the files to your computer and delete the files on the card. Then you can use the card again.

Note: The microscope supports SD card only.

3.2 Turn on the microscope

1. Press the POWER button.
2. The indicator (power LED) will light. The welcome screen will show on the LCD monitor.
3. Long press the POWER button to shutdown the microscope.

3.3 Take a Picture

1. Turn on the microscope.
2. The icon will show on the upper left corner of the screen.
3. Move the microscope to aim at the target. The target will show on the screen.
4. Hold the microscope firmly and

gently press the shutter button to take a picture.

Note:

1. User can press the ZOOM (Power) button for zooming out/in.
2. The image will be grainier as the image is digitally zoomed.

3.4 Take a Video

1. Turn on the microscope.
2. Press the FUNCTION (Right) button, select the icon. Press OK button to confirm.
3. Press the OK button again to take a video. Press the OK button again to stop.
4. Press the right button to return the function menu.

3.5 Use Digital Zoom

During taking a picture, user can press the POWER button to zoom in or zoom out. The Digital Zoom

provides 1x up to 4x magnification of your pictures.

3.6 Use Optical ZOOM

User can slide Optical Zooming switch (1) to adjust optical zoom. The optical zoom provides 3.5x and 35x magnification of the target. User can also screw the Magnification Revolving Switch to make fine adjustment.

1. Slide the switch (1) to  station, keep 60mm distance between the view case (optional) and the target. The user can view the the target in its original size.
2. Slide the switch (1) in “3.5x” station and “35x” station, the target will be magnified by 3.5x and 35x.

4. Functions

Press the Function button  to enter function menu. You can see the icons on the screen as shown. Press the Left or Right buttons to select function.

4.1 Digital video mode

Select the Digital video mode button and press the OK button to confirm. Press the OK button to take the video, and press it again to stop recording.

4.2 Digital photo mode

Select the Digital photo mode button and press the OK button to confirm. Press the OK button to take a picture.


4.3 Take Continuous Pictures

Select Take Continuous Pictures and press the OK button to confirm. The icon  will appear on the bottom left corner of the screen. Move the microscope to aim at the target. The target will show on the screen. Hold the microscope firmly and gently press the shutter button to take a picture.


Enter the function mode again, and select the Burst icon, then press the OK button to back the single picture mode.

Display Information

Digital video mode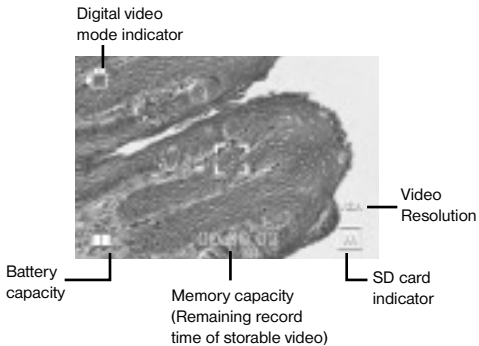

Digital photo mode

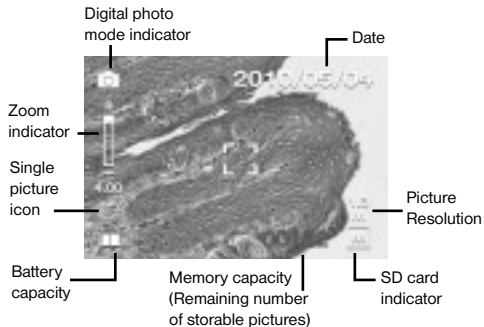4.4 Exposure value

Select EV mode and press OK button to confirm. Exposure compensation helps you increase or decrease the EV to adjust the results. This feature is applied to render the subject appropriately. The adjustment unit is 1/3, within the range -2EV ~ +2EV You can check the result of the compensation by viewing the image on the screen.

5. SETUP

In SETUP, You can see the icons shown on the screen.

5.1 Set time and date

Press the left or right button to select the date format from MM/DD/YYYY, DD/MM/YYYY or YYYY/MM/DD. Press the OK button to confirm. Then it will go to date and time setting. Press Left or Right button to adjust the date and time. Press the OK button to confirm the setting and return to main menu.

5.2 Set the picture size

Select the image quality setting to set a picture resolution. Press the left or right button to select the picture size, and press OK button to confirm and return to main menu. The Options would be 2M (1600x1200 pixels), 1.3M (1280x960 pixels), VGA (640x480 pixels).

5.3 Setting the video size

Select the video quality setting to

set the resolution. Press the left or right button to select the video size, and press OK button to confirm and return to main menu. The Options would be VGA (640x480 pixels); QVGA (320x240 pixels).

5.4 USB Connection

Select USB Connection and press OK button to confirm. PC Cam: Use your microscope as a PC camera (PC microscope). Disk Driver: Use your microscope as a mass storage for transferring files to your computer.

5.5 Light Frequency

Set the anti-flicker frequency to 50Hz. (Please use 50Hz if using the microscope in Europe, Japan and Asia). Set the anti-flicker frequency to 60 Hz. (Default setting. Please select if using the microscope in the United States).

5.6 TV Output

Connect the scanner to the TV with TV cable.

NTSC: Setting the TV Output to NTSC

PAL: Setting the TV Output to PAL

5.7 Format

Select Format and press OK button to confirm. Select '√', press OK button to format. Select 'X' and press OK button to abort.

Caution: Formatted files cannot be recovered.

5.8 Automatic Shutdown

Select Automatic Shutdown, press OK button to confirm.


5MIN: Set the microscope to power off automatically in approximately 5 minutes.

'X': Do not use this function.

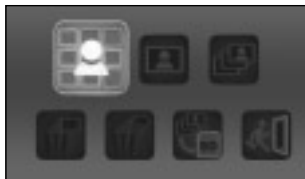
5.9 Exit

Select Exit to exit SETUP and back to Function menu.

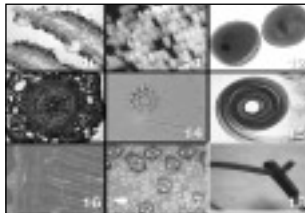6. Playback

Turn on the microscope. Press the Playback button  and you will see some pictures on the screen.

Then press the POWER button to playback mode. You will see the icons on the screen as shown. Press the Left or Right buttons to select one function.

6.1 Single

When you select the icon, press OK button to confirm. You need press Left or Right button to review the previous or next picture. Press the POWER button to go back to the Playback mode.

6.2 Slide Show

Select the Slide Show Mode and press OK button to confirm. The microscope will automatically play the photos continuously. Press any button to stop the slide show. Press the POWER button to go back to the Playback mode.

6.3 Delete One

Select the icon, press OK button and confirm with '√' to delete current image; 'X' to abort.

6.4 Delete All

Select the icon, press OK button to delete all the stored images from the internal memory or SD card. Deleted files are impossible to be recovered.

6.5 Copy to card

Select the icon, press OK button to confirm. All files will be copied into the SD card.

6.6 Exit

Select the icon, press OK button to confirm. It will return to be ready to capture image.

7. Download Files to PC

Connect the microscope to PC via USB cable. A dialogue box with options will appear as shown. You could select "Open folders to View files" to open the folder. Or select "Take no action" and find Removable disk drive in "My Computer" ("Computer" for Windows Vista) and check the folder.

8. Use the microscope as a PC-Cam

Note: Please install the microscope driver before you connect it to your PC.

Setup PC-Cam

1. Turn on the microscope and press the Function button to select SETUP. Press the OK button to confirm. Press Right / Left button to select USB Connection Icon. Select PC Cam. Press OK button to confirm.

2. Connect the microscope to PC via USB cable.

3. You could use this PC-cam with other video software you prefer. E.g. In Windows XP, Double click "my computer" (or "Computer" in Windows Vista), then click the "standard camera", you can live view in PC-Cam mode now.

9. Starting and Using the Photomizer SE Software

1. Start the Photomizer SE software on your computer.
2. For the best image result with the Photomizer SE software, we recommend setting the resolution higher. To do this, click on „adjust resolution“. If the desired resolution is not available, you can add it using the (+) field next to the resolution selection.
3. In the program window, click on „Open camera“.
4. If you have more than one device connected to your PC, you can choose the desired device in the next selection. If you only have one connected device, you do not need this step.
5. Now, you should be able to see the image from the camera on your monitor.
6. Click on „Capture“ in order to capture an image you would like to save. This will be shown on the

right in the menu bar.

7. Select this image with a single mouse click, then click on „Transfer image“.
8. You will now leave the image-capture stage, and go back to the program window.
9. Now click at the top of the screen on “File” and then on “save file” in order to secure the image.

Tip:

If you have questions or problems with the “Photomizer SE” software, you can always go to the program’s menu bar, and click on „?“ and then on „Open help“. For questions or problems, you can look on the website www.photomizer.net. Please ask your parents before using the Internet.

Technical Data	
Sensor Resolution	8.5 mm (1 / 3") CMOS, 1280 x 1024 Pixel
Lens Specification	F8.0, f = 7.1 mm, View Angle 44°, 2G2P
Illumination	Built-in LED x 4
O / S	Microsoft Windows XP/ Vista / 7
Magnification	6 cm (2.4") Screen: 3.5x / 35x 43 cm (17") Screen: 19x / 190x
White Balance	Auto
Exposure Control	Automatic
Capacity Still Mode	VGA: 640 x 480 Pixel 1.3M: 1280 x 960 Pixel 2M: 1600 x 1200 Pixel
TV OUT	PAL / NTSC (Optional)
Video Mode	Video: (w/o Audio function) VGA 640 x 480 Pixel, 30 fps QVGA 320 x 240 Pixel, 19 fps
Compression	JPEG / MJPEG (AVI)
Digital zoom	1x - 4 x Digital Zoom for Preview & Playback
Display	6cm (2.4") LCD MONITOR

PC Interface	USB 1.1, MSD Funktion
Storage Type	SD Card Slot (SD card not included)
Power	2 x "AA" 1.5V, Alkaline Batteries
Dimensions (LxWxH)	54.5x73x96.5 mm 40x64.5x15 mm (view case)

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